

ENG

Instructions for Use:
HUMAN sCD40 ELISA

Catalogue number:
RAF092R

For research use only!

 **BioVendor**
R&D[®]



BioVendor – Laboratorní medicína a.s.

Karásek 1767/1, 621 00 Brno, Czech Republic

+420 549 124 185

info@biovendor.com

sales@biovendor.com

www.biovendor.com

1. INTENDED USE	3
2. SUMMARY	3
3. PRINCIPLES OF THE TEST	4
4. REAGENTS PROVIDED	5
5. STORAGE INSTRUCTIONS – ELISA KIT	5
6. SPECIMEN COLLECTION AND STORAGE INSTRUCTIONS	6
7. MATERIALS REQUIRED BUT NOT PROVIDED	6
8. PRECAUTIONS FOR USE	6
9. PREPARATION OF REAGENTS	7
10. TEST PROTOCOL	9
11. CALCULATION OF RESULTS	11
12. LIMITATIONS	14
13. PERFORMANCE CHARACTERISTICS	14
14. REAGENT PREPARATION SUMMARY	17
15. TEST PROTOCOL SUMMARY	18
16. EXPLANATION OF SYMBOLS	19

HISTORY OF CHANGES

Previous version	Current version
ENG.004.A	ENG.005.A
Formal change of document appearance (rebranding).	

1. INTENDED USE

The Human sCD40 ELISA is an enzyme-linked immunosorbent assay for the quantitative detection of Human CD40.

2. SUMMARY


CD40 is a 50 kDa membrane-bound type I glycoprotein expressed by numerous cells, most notably B lymphocytes and monocytes and antigen presenting cells (APC) such as macrophages, dendritic cells and fibroblasts. It is furthermore highly expressed on various malignant cells. The expression of CD40 regulates T-cell - APC interaction and has been shown to be centrally involved in a wide array of inflammatory events. The function of CD40 is very broad. CD40 belongs to the TNF-receptor family. The ligand for CD40 (CD154) is a 33 kDa type II transmembrane protein mainly expressed by activated T-cells, and is a member of the TNF superfamily. CD40/CD40L interactions are essential for T-cell-dependent B cell proliferation and differentiation, for activation of antigen presenting cells and for cytokine production on numerous other cells.

It is presently accepted that CD40 plays a critical role in the regulation of immune responses.

3. PRINCIPLES OF THE TEST


An anti-Human CD40 coating antibody is adsorbed onto microwells.

Figure 1 Coated Microwell

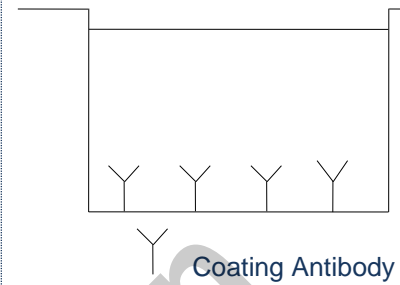


Human CD40 present in the sample or standard binds to antibodies adsorbed to the microwells. A biotin-conjugated anti-Human CD40 antibody is added and binds to Human CD40 captured by the first antibody.

Figure 2 First Incubation

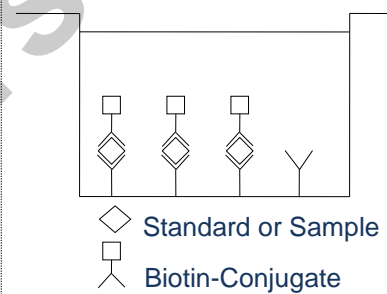


Following incubation unbound biotin-conjugated anti-Human CD40 antibody is removed during a wash step. Streptavidin-HRP is added and binds to the biotin-conjugated anti-Human CD40 antibody.

Figure 3 Second Incubation

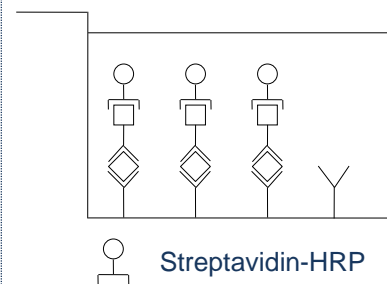


Following incubation unbound Streptavidin-HRP is removed during a wash step, and substrate solution reactive with HRP is added to the wells.

Figure 4 Third Incubation

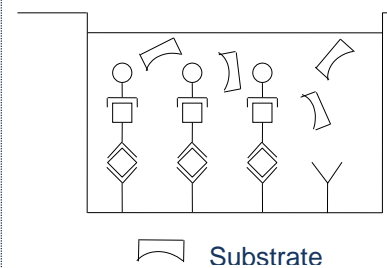

A coloured product is formed in proportion to the amount of Human CD40 present in the sample or standard. The reaction is terminated by addition of acid and absorbance is measured at 450 nm. A standard curve is prepared from 7 Human CD40 standard dilutions and Human CD40 sample concentration determined.

Figure 5

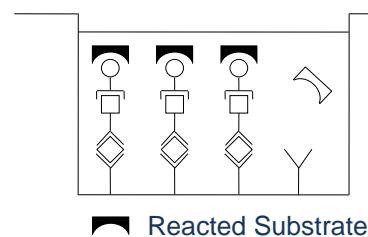

4. REAGENTS PROVIDED

- 1 aluminium pouch with a Antibody Coated Microtiter Strips with monoclonal antibody to Human CD40
- 1 vial (0,07 ml) Biotin-Conjugate anti-Human CD40 monoclonal antibody
- 1 vial (0,15 ml) Streptavidin-HRP
- 2 vials Human CD40 Standard lyophilized, 1 ng/ml upon reconstitution
- 1 vial (12 ml) Sample Diluent
- 1 vial (5 ml) Assay Buffer Concentrate 20x (PBS with 1% Tween 20 and 10% BSA)
- 1 bottle (50 ml) Wash Buffer Concentrate 20x (PBS with 1% Tween 20)
- 1 vial (15 ml) Substrate Solution (tetramethyl-benzidine)
- 1 vial (15 ml) Stop Solution (1M Phosphoric acid)
- 4 Adhesive Films

5. STORAGE INSTRUCTIONS – ELISA KIT

Store kit reagents between 2° and 8°C. Immediately after use remaining reagents should be returned to cold storage (2° to 8°C). Expiry of the kit and reagents is stated on labels.

Expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, this reagent is not contaminated by the first handling.

6. SPECIMEN COLLECTION AND STORAGE INSTRUCTIONS

Cell culture supernatant, serum, plasma (EDTA, heparin) and amniotic fluid were tested with this assay. Other biological samples might be suitable for use in the assay. Remove serum or plasma from the clot or cells as soon as possible after clotting and separation.

Samples containing a visible precipitate must be clarified prior to use in the assay. Do not use grossly hemolyzed or lipemic specimens.

Samples should be aliquoted and must be stored frozen at -20°C to avoid loss of bioactive Human CD40. If samples are to be run within 24 hours, they may be stored at 2° to 8°C (for sample stability refer to 13.5).

Avoid repeated freeze-thaw cycles. Prior to assay, the frozen sample should be brought to room temperature slowly and mixed gently.

7. MATERIALS REQUIRED BUT NOT PROVIDED

- 5 ml and 10 ml graduated pipettes
- 5 μl to 1000 μl adjustable single channel micropipettes with disposable tips
- 50 μl to 300 μl adjustable multichannel micropipette with disposable tips
- Multichannel micropipette reservoir
- Beakers, flasks, cylinders necessary for preparation of reagents
- Device for delivery of wash solution (multichannel wash bottle or automatic wash system)
- Microwell strip reader capable of reading at 450 nm (620 nm as optional reference wave length)
- Glass-distilled or deionized water
- Statistical calculator with program to perform regression analysis

8. PRECAUTIONS FOR USE

- All chemicals should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statement(s) for specific advice.
- Reagents are intended for research use only and are not for use in diagnostic or therapeutic procedures.
- Do not mix or substitute reagents with those from other lots or other sources.
- Do not use kit reagents beyond expiration date on label.
- Do not expose kit reagents to strong light during storage or incubation.
- Do not pipette by mouth.
- Do not eat or smoke in areas where kit reagents or samples are handled.
- Avoid contact of skin or mucous membranes with kit reagents or specimens.
- Rubber or disposable latex gloves should be worn while handling kit reagents or specimens.
- Avoid contact of substrate solution with oxidizing agents and metal.
- Avoid splashing or generation of aerosols.
- In order to avoid microbial contamination or cross-contamination of reagents or specimens which may invalidate the test use disposable pipette tips and/or pipettes.
- Use clean, dedicated reagent trays for dispensing the conjugate and substrate reagent.

- Exposure to acid inactivates the conjugate.
- Glass-distilled water or deionized water must be used for reagent preparation.
- Substrate solution must be at room temperature prior to use.
- Decontaminate and dispose specimens and all potentially contaminated materials as they could contain infectious agents. The preferred method of decontamination is autoclaving for a minimum of 1 hour at 121.5°C.
- Liquid wastes not containing acid and neutralized waste may be mixed with sodium hypochlorite in volumes such that the final mixture contains 1.0% sodium hypochlorite. Allow 30 minutes for effective decontamination. Liquid waste containing acid must be neutralized prior to the addition of sodium hypochlorite.

9. PREPARATION OF REAGENTS

Buffer Concentrates should be brought to room temperature and should be diluted before starting the test procedure.

If crystals have formed in the **Buffer Concentrates**, warm them gently until they have completely dissolved.

9.1 Wash Buffer (1x)

Pour entire contents (50 ml) of the **Wash Buffer Concentrate** (20x) into a clean 1000 ml graduated cylinder. Bring to final volume of 1000 ml with glass-distilled or deionized water. Mix gently to avoid foaming.

Transfer to a clean wash bottle and store at 2° to 25°C. Please note that Wash Buffer (1x) is stable for 30 days.

Wash Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Wash Buffer Concentrate (20x) (ml)	Distilled Water (ml)
1 - 6	25	475
1 - 12	50	950

9.2 Assay Buffer (1x)

Pour the entire contents (5 ml) of the **Assay Buffer Concentrate** (20x) into a clean 100 ml graduated cylinder. Bring to final volume of 100 ml with distilled water. Mix gently to avoid foaming.

Store at 2° to 8°C. Please note that the Assay Buffer (1x) is stable for 30 days.

Assay Buffer (1x) may also be prepared as needed according to the following table:

Number of Strips	Assay Buffer Concentrate (20x) (ml)	Distilled Water (ml)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

9.3 Biotin-Conjugate

Please note that the **Biotin-Conjugate** should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated Biotin-Conjugate solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Biotin-Conjugate (ml)	Assay Buffer (1x) (ml)
1 - 6	0.03	2.97
1 - 12	0.06	5.94

9.4 Streptavidin-HRP

Please note that the **Streptavidin-HRP** should be used within 30 minutes after dilution.

Make a 1:100 dilution of the concentrated Streptavidin-HRP solution with Assay Buffer (1x) in a clean plastic tube as needed according to the following table:

Number of Strips	Streptavidin-HRP (ml)	Assay Buffer (1x) (ml)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

9.5 Human CD40 Standard

Reconstitute **Human CD40 standard** by addition distilled water. Reconstitution volume is stated in the Quality Control Sheet. Swirl or mix gently to insure complete and homogeneous solubilization (concentration of reconstituted standard = 1000 pg/ml).

Allow the reconstituted standard to sit for 10-30 minutes. Mix well prior to making dilutions.

After usage remaining standard cannot be stored and has to be discarded.

9.5.1 External Standard Dilution

Label 7 tubes, one for each standard point.

S1, S2, S3, S4, S5, S6, S7


Then prepare 1:2 serial dilutions for the standard curve as follows:

Pipette 225 µl of Sample Diluent into each tube.

Pipette 225 µl of reconstituted standard (concentration = 1000 pg/ml) into the first tube, labelled S1, and mix (concentration of standard 1 = 500 pg/ml). Pipette 225 µl of this dilution into the second tube, labelled S2, and mix thoroughly before the next transfer. Repeat serial dilutions 5 more times thus creating the points of the standard curve (see Figure 6).

Sample Diluent serves as blank.


Figure 6

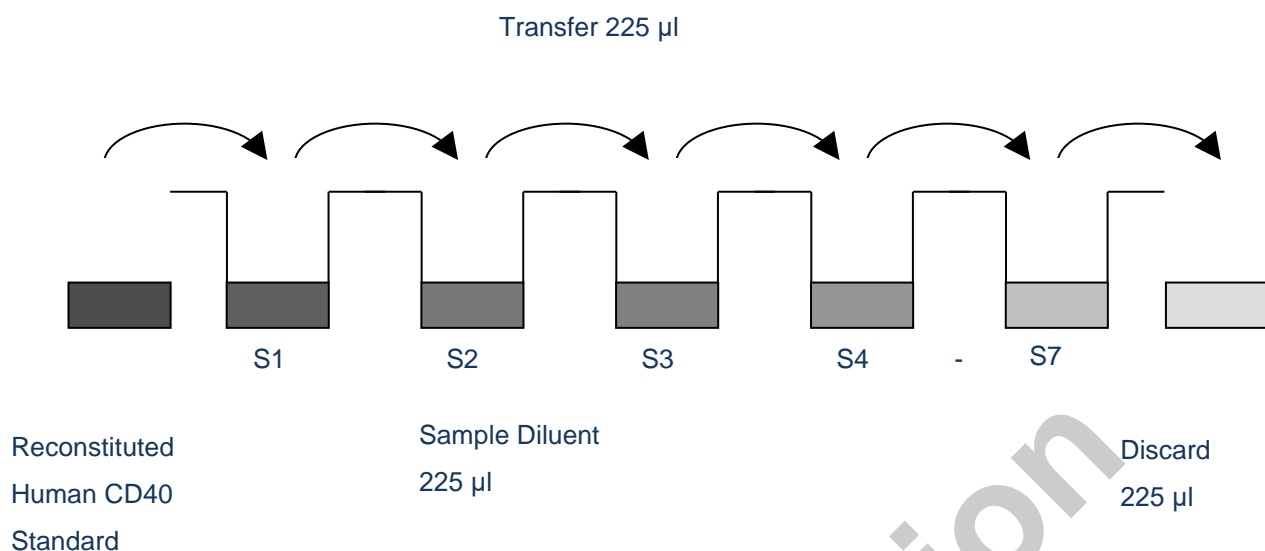

10. TEST PROTOCOL

- Determine the number of microwell strips required to test the desired number of samples plus appropriate number of wells needed for running blanks and standards. Each sample, standard, blank and optional control sample should be assayed in duplicate. Remove extra microwell strips from holder and store in foil bag with the desiccant provided at 2°-8°C sealed tightly.
- Wash the microwell strips twice with approximately 400 μ l **Wash Buffer** per well with thorough aspiration of microwell contents between washes. Allow the Wash Buffer to sit in the wells for about **10 – 15 seconds** before aspiration. Take care not to scratch the surface of the microwells. After the last wash step, empty wells and tap microwell strips on absorbent pad or paper towel to remove excess Wash Buffer. Use the microwell strips immediately after washing. Alternatively microwell strips can be placed upside down on a wet absorbent paper for not longer than 15 minutes. **Do not allow wells to dry.**
- Standard dilution on the microwell plate** (Alternatively the standard dilution can be prepared in tubes – see 9.5.1): Add 100 μ l of Sample Diluent in duplicate to all **standard wells**. Pipette 100 μ l of prepared **standard** (see Preparation of Standard 9.5., concentration = 1000 pg/ml) in duplicate into well A1 and A2 (see Table 1). Mix the contents of wells A1 and A2 by repeated aspiration and ejection (concentration of standard 1, S1 = 500 pg/ml), and transfer 100 μ l to wells B1 and B2, respectively (see Figure 7). Take care not to scratch the inner surface of the microwells. Continue this procedure 5 times, creating two rows of Human CD40 standard dilutions ranging from 500.0 to 7.8 pg/ml. Discard 100 μ l of the contents from the last microwells (G1, G2) used.

Figure 7

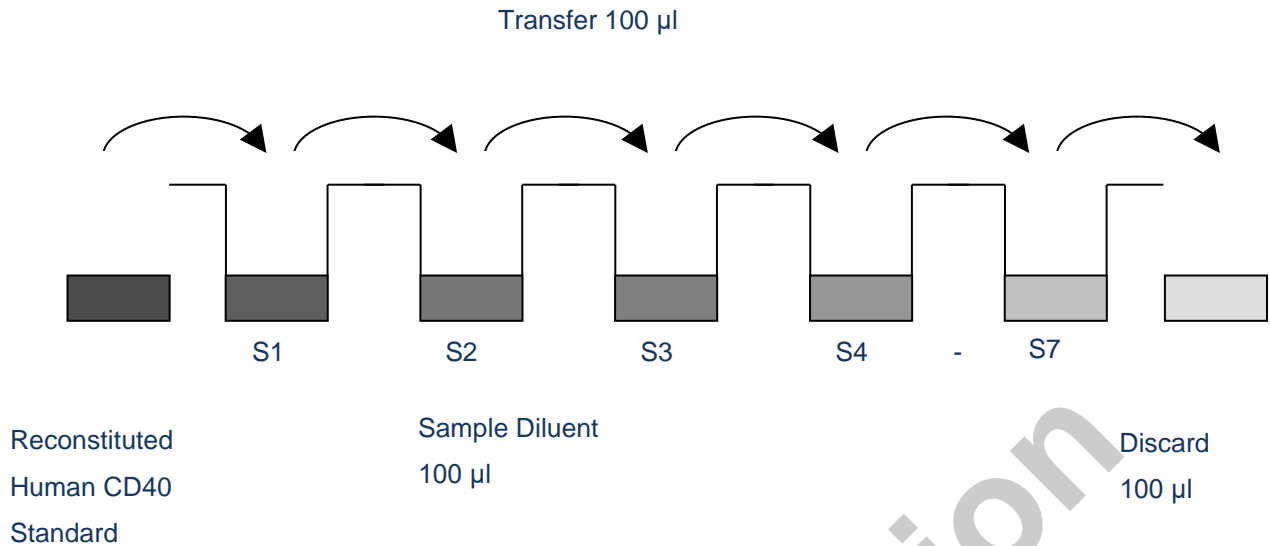

In case of an **external standard dilution**, pipette 100 μ l of these standard dilutions (S1 - S7) in the standard wells according to Table 1.

Table 1 Table depicting an example of the arrangement of blanks, standards and samples in the microwell strips:

	1	2	3	4
A	Standard 1 (500.0 pg/ml)	Standard 1 (500.0 pg/ml)	Sample 1	Sample 1
B	Standard 2 (250.0 pg/ml)	Standard 2 (250.0 pg/ml)	Sample 2	Sample 2
C	Standard 3 (125.0 pg/ml)	Standard 3 (125.0 pg/ml)	Sample 3	Sample 3
D	Standard 4 (62.5 pg/ml)	Standard 4 (62.5 pg/ml)	Sample 4	Sample 4
E	Standard 5 (31.3 pg/ml)	Standard 5 (31.3 pg/ml)	Sample 5	Sample 5
F	Standard 6 (15.6 pg/ml)	Standard 6 (15.6 pg/ml)	Sample 6	Sample 6
G	Standard 7 (7.8 pg/ml)	Standard 7 (7.8 pg/ml)	Sample 7	Sample 7
H	Blank	Blank	Sample 8	Sample 8

4. Add 100 μ l of **Sample Diluent** in duplicate to the **blank wells**.
5. Add 50 μ l of **Sample Diluent** to the **sample wells**.
6. Add 50 μ l of each **sample** in duplicate to the **sample wells**.
7. Prepare **Biotin-Conjugate** (see Preparation of Biotin-Conjugate 9.3.).
8. Add 50 μ l of **Biotin-Conjugate** to all wells.
9. Cover with an adhesive film and incubate at room temperature (18 to 25°C) for 2 hours, if available on a microplate shaker.
10. Prepare **Streptavidin-HRP** (refer to Preparation of Streptavidin-HRP 9.4.).
11. Remove adhesive film and empty wells. **Wash** microwell strips 6 times according to point 2. of the test protocol. Proceed immediately to the next step.

12. Add 100 µl of diluted **Streptavidin-HRP** to all wells, including the blank wells.
13. Cover with an adhesive film and incubate at room temperature (18° to 25°C) for 1 hour, if available on a microplate shaker.
14. Remove adhesive film and empty wells. **Wash** microwell strips 6 times according to point 2. of the test protocol. Proceed immediately to the next step.
15. Pipette 100 µl of **TMB Substrate Solution** to all wells.
16. Incubate the microwell strips at room temperature (18° to 25°C) for about 10 min. Avoid direct exposure to intense light. **The colour development on the plate should be monitored and the substrate reaction stopped (see next point of this protocol) before positive wells are no longer properly recordable. Determination of the ideal time period for colour development has to be done individually for each assay.** It is recommended to add the stop solution when the highest standard has developed a dark blue colour. Alternatively the colour development can be monitored by the ELISA reader at 620 nm. The substrate reaction should be stopped as soon as Standard 1 has reached an OD of 0.9 – 0.95.
17. Stop the enzyme reaction by quickly pipetting 100 µl of **Stop Solution** into each well. It is important that the Stop Solution is spread quickly and uniformly throughout the microwells to completely inactivate the enzyme. Results must be read immediately after the Stop Solution is added or within one hour if the microwell strips are stored at 2 - 8°C in the dark.
18. Read absorbance of each microwell on a spectro-photometer using 450 nm as the primary wave length (optionally 620 nm as the reference wave length; 610 nm to 650 nm is acceptable). Blank the plate reader according to the manufacturer's instructions by using the blank wells. Determine the absorbance of both the samples and the standards.

Note: In case of incubation without shaking the obtained O.D. values may be lower than indicated below. Nevertheless the results are still valid.

11. CALCULATION OF RESULTS

- Calculate the average absorbance values for each set of duplicate standards and samples. Duplicates should be within 20 per cent of the mean value.
- Create a standard curve by plotting the mean absorbance for each standard concentration on the ordinate against the Human CD40 concentration on the abscissa. Draw a best fit curve through the points of the graph (a 5-parameter curve fit is recommended).
- To determine the concentration of circulating Human CD40 for each sample, first find the mean absorbance value on the ordinate and extend a horizontal line to the standard curve. At the point of intersection, extend a vertical line to the abscissa and read the corresponding Human CD40 concentration.
- If instructions in this protocol have been followed samples have been diluted 1:2 (50 µl sample + 50 µl Sample Diluent), the concentration read from the standard curve must be multiplied by the dilution factor (x 2).
- Calculation of samples with a concentration exceeding standard 1 may result in incorrect, low Human CD40 levels. Such samples require further external predilution according to expected Human CD40 values with Sample Diluent in order to precisely quantitate the actual Human CD40 level.
- It is suggested that each testing facility establishes a control sample of known Human CD40 concentration and runs this additional control with each assay. If the values obtained are not within the expected range of the control, the assay results may be invalid.
- A representative standard curve is shown in Figure 8. This curve cannot be used to derive test results. Each laboratory must prepare a standard curve for each group of microwell strips assayed.

Figure 8 Representative standard curve for Human CD40 ELISA. Human CD40 was diluted in serial 2-fold steps in Sample Diluent.

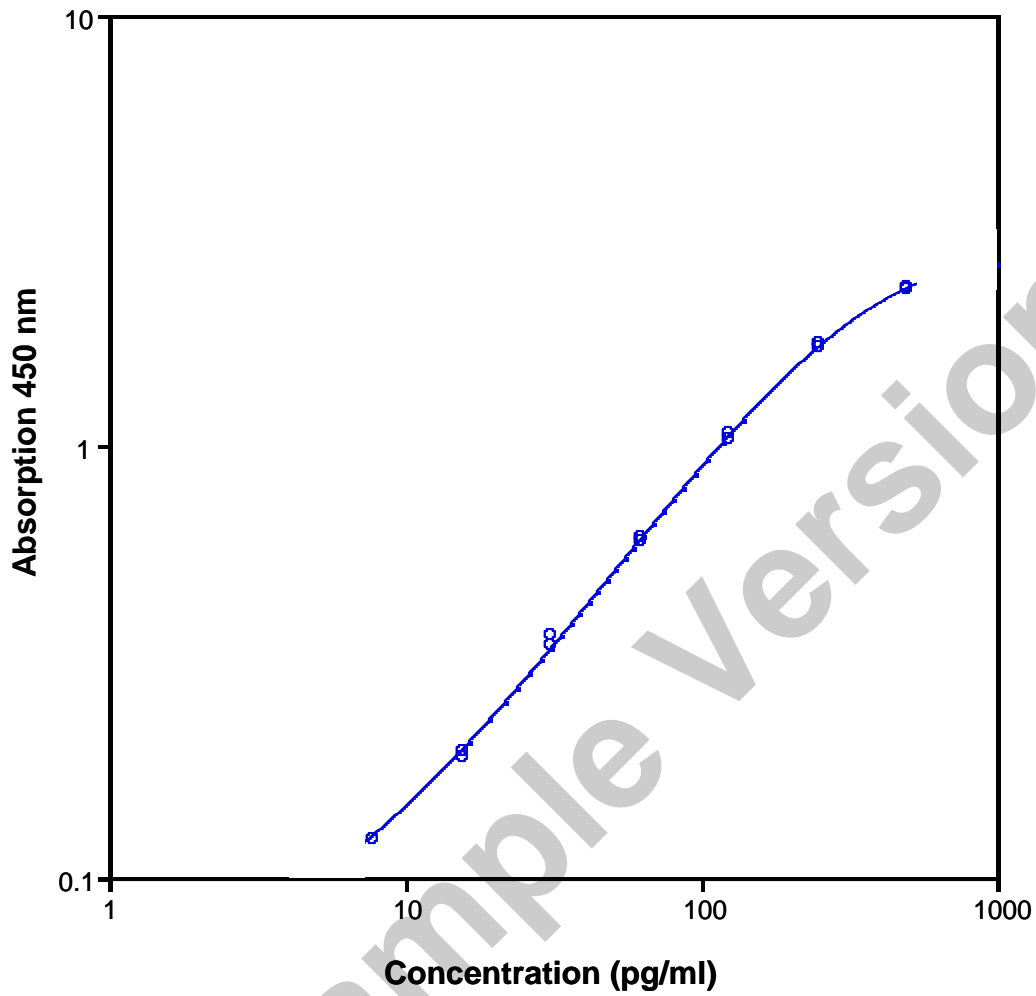


Table 2 Typical data using the Human sCD40 ELISA

Measuring wavelength: 450 nm
 Reference wavelength: 620 nm

Standard	Human sCD40 Concentration (pg/ml)	O.D. at 450 nm	Mean O.D. at 450 nm	C.V.(%)
1	500.0	2.325	2.330	0.3
		2.334		
2	250.0	1.727	1.711	1.4
		1.694		
3	125.0	1.069	1.052	2.4
		1.034		
4	62.5	0.604	0.610	1.4
		0.616		
5	31.3	0.345	0.354	3.6
		0.363		
6	15.6	0.197	0.193	2.9
		0.189		
7	7.8	0.122	0.122	0.0
		0.122		
Blank	0.0	0.024	0.028	20.2
		0.032		

The OD values of the standard curve may vary according to the conditions of assay performance (e.g. operator, pipetting technique, washing technique or temperature effects). Furthermore shelf life of the kit may affect enzymatic activity and thus colour intensity. Values measured are still valid.

12. LIMITATIONS

- Since exact conditions may vary from assay to assay, a standard curve must be established for every run.
- Bacterial or fungal contamination of either screen samples or reagents or cross-contamination between reagents may cause erroneous results.
- Disposable pipette tips, flasks or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.
- Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Empty wells completely before dispensing fresh wash solution, fill with Wash Buffer as indicated for each wash cycle and do not allow wells to sit uncovered or dry for extended periods.
- The use of radioimmunotherapy has significantly increased the number of patients with Human anti-mouse IgG antibodies (HAMA). HAMA may interfere with assays utilizing murine monoclonal antibodies leading to both false positive and false negative results. Serum samples containing antibodies to murine immunoglobulins can still be analysed in such assays when murine immunoglobulins (serum, ascitic fluid, or monoclonal antibodies of irrelevant specificity) are added to the sample.

13. PERFORMANCE CHARACTERISTICS

13.1 Sensitivity

The limit of detection of Human CD40 defined as the analyte concentration resulting in an absorbance significantly higher than that of the dilution medium (mean plus 2 standard deviations) was determined to be 1.3 pg/ml (mean of 6 independent assays).

13.2 Reproducibility

13.2.1 Intra-assay

Reproducibility within the assay was evaluated in 2 independent experiments. Each assay was carried out with 4 replicates of 8 serum samples containing different concentrations of Human CD40. 2 standard curves were run on each plate. Data below show the mean Human CD40 concentration and the coefficient of variation for each sample (see Table 3). The calculated overall intra-assay coefficient of variation was 5.5%.

Table 3 The mean Human CD40 concentration and the coefficient of variation for each sample

Sample	Experiment	Mean Human CD40 Concentration (pg/ml)	Coefficient of Variation (%)
1	1	221.2	8.7
	2	224.9	2.0
2	1	105.2	3.0
	2	133.8	8.7
3	1	272.8	4.6
	2	269.6	0.7
4	1	223.2	0.5
	2	222.8	7.9

Sample	Experiment	Mean Human CD40 Concentration (pg/ml)	Coefficient of Variation (%)
5	1	51.6	11.1
	2	58.3	12.5
6	1	153.3	13.0
	2	138.8	8.3
7	1	1630.1	3.3
	2	1308.8	0.4
8	1	742.3	0.4
	2	682.5	2.9

13.2.2 Inter-assay

Assay to assay reproducibility within one laboratory was evaluated in 2 independent experiments. Each assay was carried out with 4 replicates of 8 serum samples containing different concentrations of Human CD40. 2 standard curves were run on each plate. Data below show the mean Human CD40 concentration and the coefficient of variation calculated on 8 determinations of each sample (see Table 4). The calculated overall inter-assay coefficient of variation was 7.0%.

Table 4 The mean Human CD40 concentration and the coefficient of variation of each sample

Sample	Mean Human sCD40 Concentration (pg/ml)	Coefficient of Variation (%)
1	223.0	1.2
2	119.5	16.9
3	271.2	0.8
4	223.0	0.1
5	54.9	8.7
6	149.1	7.0
7	1469.4	15.5
8	712.4	5.9

13.3 Spiking Recovery

The spiking recovery was evaluated by spiking 4 levels of Human CD40 into pooled normal serum samples. Recoveries were determined in 3 independent experiments with 6 replicates each. The amount of endogenous Human CD40 in unspiked serum was subtracted from the spike values. The recovery ranged from 81% to 110% with an overall mean recovery of 94%.

13.4 Dilution Linearity

4 serum samples with different levels of Human CD40 were analysed at serial 2 fold dilutions with 4 replicates each.

The recovery ranged from 82% to 104% with an overall recovery of 94% (see Table 5).

Table 5

Sample	Dilution	Expected Human CD40 Concentration (pg/ml)	Observed Human CD40 Concentration (pg/ml)	Recovery of Expected Human CD40 Concentration (%)
1	1:2	-	121.4	-
	1:4	60.7	62.9	104
	1:8	30.3	31.4	104
2	1:2	-	88.0	-
	1:4	44.0	43.2	98
	1:8	22.0	21.4	97
3	1:2	-	419.9	-
	1:4	209.9	197.2	94
	1:8	105.0	85.7	82
4	1:2	-	282.0	-
	1:4	141.0	130.7	93
	1:8	70.5	58.8	83

13.5 Sample Stability

13.5.1 Freeze-Thaw Stability

Aliquots of serum samples (spiked or unspiked) were stored at -20°C and thawed 5 times, and the Human CD40 levels determined. There was no significant loss of Human CD40 immunoreactivity detected by freezing and thawing.

13.5.2 Storage Stability

Aliquots of serum samples (spiked or unspiked) were stored at -20°C, 2-8°C, room temperature (RT) and at 37°C, and the Human CD40 level determined after 24 h. There was no significant loss of Human sCD40 immunoreactivity detected during storage under above conditions.

13.6 Comparison of Serum and Plasma

From 8 individuals, serum as well as EDTA, citrate and heparin plasma obtained at the same time point were evaluated. Human CD40 concentrations were not significantly different and therefore all these body fluids are suitable for the assay. It is nevertheless highly recommended to assure the uniformity of blood preparations.

13.7 Specificity

The interference of circulating factors of the immune system was evaluated by spiking these proteins at physiologically relevant concentrations into a Human CD40 positive serum. There was no crossreactivity detected.

13.8 Expected Values

A panel of 8 serum samples from randomly selected apparently healthy donors was tested for Human CD40.

The detected Human CD40 levels ranged between 53.3 and 156.9 pg/ml with a mean level of 102.1 pg/ml and a standard deviation of 37.1 pg/ml.

14. REAGENT PREPARATION SUMMARY

14.1 Wash Buffer (1x)

Add **Wash Buffer Concentrate** 20x (50 ml) to 950 ml distilled water.

Number of Strips	Wash Buffer Concentrate (ml)	Distilled Water (ml)
1 - 6	25	475
1 - 12	50	950

14.2 Assay Buffer (1x)

Add **Assay Buffer Concentrate** 20x (5 ml) to 95 ml distilled water.

Number of Strips	Assay Buffer Concentrate (ml)	Distilled Water (ml)
1 - 6	2.5	47.5
1 - 12	5.0	95.0

14.3 Biotin-Conjugate

Make a 1:100 dilution of **Biotin-Conjugate** in Assay Buffer (1x):

Number of Strips	Biotin-Conjugate (ml)	Assay Buffer (1x) (ml)
1 - 6	0.03	2.97
1 - 12	0.06	5.94

14.4 Streptavidin-HRP

Make a 1:100 dilution of **Streptavidin-HRP** in Assay Buffer (1x):

Number of Strips	Streptavidin-HRP (ml)	Assay Buffer (1x) (ml)
1 - 6	0.06	5.94
1 - 12	0.12	11.88

14.5 Human CD40 Standard










Reconstitute lyophilized **Human CD40 standard** with distilled water. (Reconstitution volume is stated in the Quality Control Sheet).

15. TEST PROTOCOL SUMMARY

1. Determine the number of microwell strips required.
2. Wash microwell strips twice with Wash Buffer.
3. Standard dilution on the microwell plate: Add 100 µl Sample Diluent, in duplicate, to all standard wells. Pipette 100 µl prepared standard into the first wells and create standard dilutions by transferring 100 µl from well to well. Discard 100 µl from the last wells. Alternatively external standard dilution in tubes (see 9.5.1): Pipette 100 µl of these standard dilutions in the microwell strips.
4. Add 100 µl Sample Diluent, in duplicate, to the blank wells.
5. Add 50 µl Sample Diluent to sample wells.
6. Add 50 µl sample in duplicate, to designated sample wells.
7. Prepare Biotin-Conjugate.
8. Add 50 µl Biotin-Conjugate to all wells.
9. Cover microwell strips and incubate 2 hours at room temperature (18° to 25°C).
10. Prepare Streptavidin-HRP.
11. Empty and wash microwell strips 6 times with Wash Buffer.
12. Add 100 µl diluted Streptavidin-HRP to all wells.
13. Cover microwell strips and incubate 1 hour at room temperature (18° to 25°C).
14. Empty and wash microwell strips 6 times with Wash Buffer.
15. Add 100 µl of TMB Substrate Solution to all wells.
16. Incubate the microwell strips for about 10 minutes at room temperature (18° to 25°C).
17. Add 100 µl Stop Solution to all wells.
18. Blank microwell reader and measure colour intensity at 450 nm.

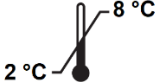Note: If instructions in this protocol have been followed samples have been diluted 1:2 (50 µl sample + 50 µl Sample Diluent), the concentration read from the standard curve must be multiplied by the dilution factor (x 2).

16. EXPLANATION OF SYMBOLS

	Catalogue number
	Batch code
	Caution
	Use by date
	Temperature limit
	Manufacturer
 <p data-bbox="260 1184 467 1216">www.biovendor.com</p>	Read electronic instructions for use - eIFU
	The content is sufficient for 96 tests
	Biological risks



BioVendor – Laboratorní medicína a.s.

Karásek 1767/1, 621 00 Brno, Czech Republic

+420 549 124 185

info@biovendor.com

sales@biovendor.com

www.biovendor.com



FOLLOW US