


MicroRNA-BASED DIAGNOSTICS IN CARDIOVASCULAR DISEASE

J. Jancovicova¹, M. Hlozankova¹, B. Dvorakova¹, M. Buresova¹, M. Hamanova¹, L. Chalupova¹,
K. Cuchnova¹, A. Kappel², A. Keller³, V. Ruzicka¹

¹ BioVendor-Laboratorní medicína a.s., Research & Diagnostic Division, Brno, Czech Republic

² Corporate Technology, Siemens AG, Erlangen, Germany

³ Clinical Bioinformatics, Medical Faculty, Saarland University, Germany

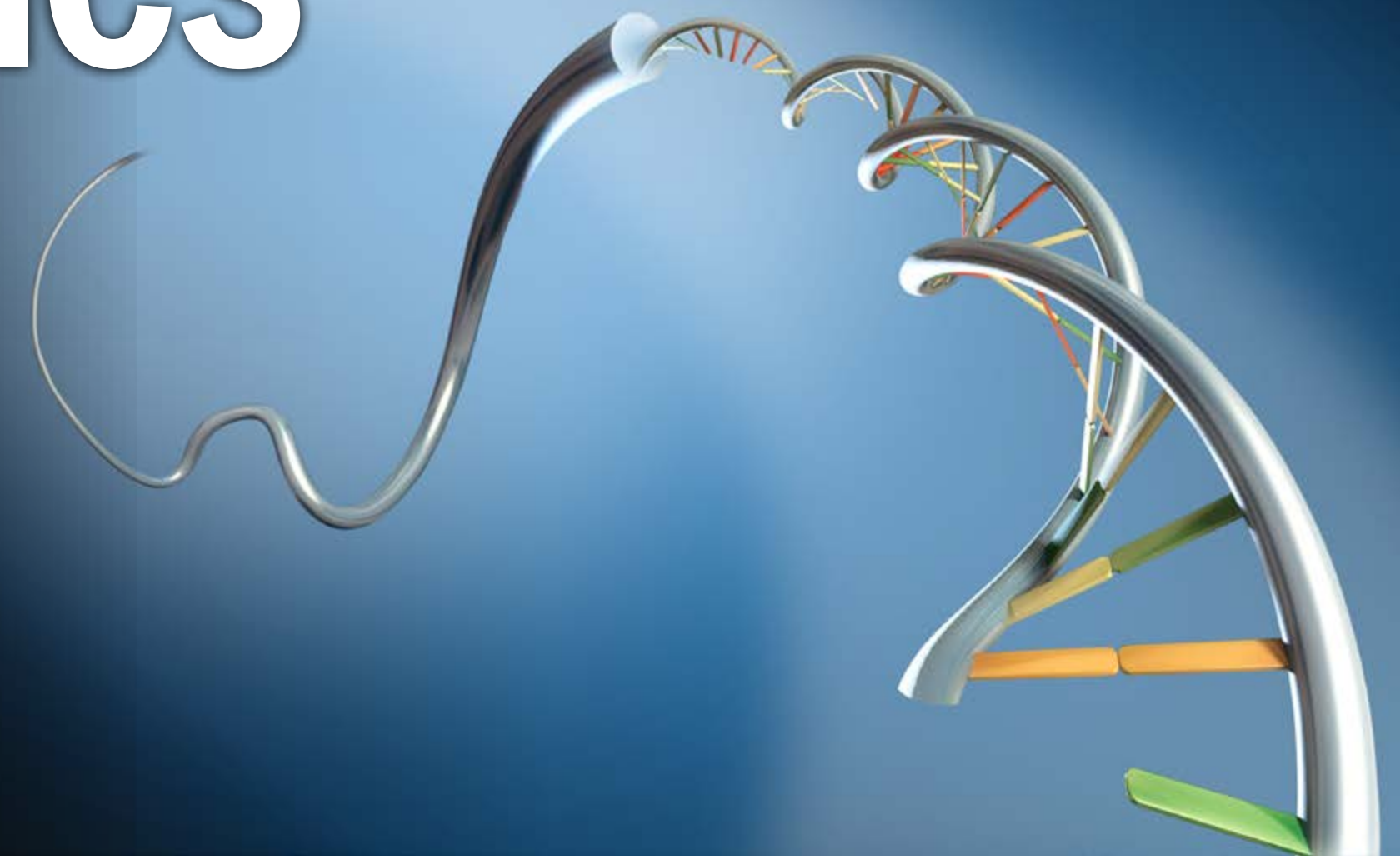

Introduction

MicroRNAs (miRNAs) are small non-coding RNA molecules playing an important regulatory role in gene translation through silencing or degradation of target mRNAs. They are involved in a wide range of biological processes, including differentiation and proliferation, metabolism, hemostasis, apoptosis or inflammation, and also in pathophysiology of many diseases. Recently, numerous studies have suggested circulating miRNAs as promising diagnostic and prognostic biomarkers of many diseases.

MicroRNA-21 (miR-21) is highly expressed in the cardiovascular system. Recent studies have revealed that its expression is dysregulated in heart under cardiovascular disease conditions such as proliferative vascular disease, cardiac hypertrophy and heart failure or ischemic heart disease. miR-21 has been found to play important roles in vascular smooth muscle cell proliferation and apoptosis, cardiac cell growth and death and cardiac fibroblast functions.

Method


We are introducing a novel, immunoassay-based method of miRNA quantification which involves hybridization of miRNA isolated from a patient sample to complementary biotinylated DNA oligonucleotide probe. The DNA/RNA hybrids are then transferred onto a stationary solid phase coated with monoclonal antibody specific to perfectly matched DNA/miRNA hybrids. After washing, the solid phase is incubated with streptavidin-HRP conjugate and the resulting complexes are visualized (after another washing step) by a chromogenic substrate. Our immunoassay exhibits superior analytical specificity, limit of detection as low as 0.1 attomol/μl miRNA, excellent analytical characteristics and strong correlation with the qRT-PCR method (Pearson correlation coefficient >0.9).

Moreover, the assay can be run on common immunoassay analyzers, is compatible with standard clinical workflow, does not require amplification steps and results are obtained in less than three hours including miRNA profiling. Our method enables to analyze miRNA using conventional immunoassay analyzers, thus, it can promote utilization of miRNA biomarkers in clinical and laboratory practice.

References:


- Shenoy, Archana, and Robert H. Bleiloch. *Regulation of microRNA function in somatic stem cell proliferation and differentiation*. Nature Reviews Molecular Cell Biology 15.9 (2014): 565-576.
- Rottiers, Veerle, and Anders M. Näär. *MicroRNAs in metabolism and metabolic disorders*. Nature reviews Molecular cell biology 13.4 (2012): 239-250.
- Teruel-Montoya, R., F. R. Rosendaal, and C. Martínez. *MicroRNAs in hemostasis*. Journal of Thrombosis and Haemostasis 13.2 (2015): 170-181.
- Jovanovic, M., and M. O. Hengartner. *miRNAs and apoptosis: RNAs to die for*. Oncogene 25.46 (2006): 6176-6187.
- Singh, Ram Pyare, et al. *The role of miRNA in inflammation and autoimmunity*. Autoimmunity reviews 12.12 (2013): 1160-1165.
- Esquele-Kerscher, Aurora, and Frank J. Slack. *Oncomirs—microRNAs with a role in cancer*. Nature Reviews Cancer 6.4 (2006): 259-269.
- Romaine, Simon PR, et al. *MicroRNAs in cardiovascular disease: an introduction for clinicians*. Heart 101.12 (2015): 921-928.
- Tan, Lin, et al. *Genome-wide serum microRNA expression profiling identifies serum biomarkers for Alzheimer's disease*. Journal of Alzheimer's Disease 40.4 (2014): 1017-1027.
- Wang, Jin, Jinyun Chen, and Subrata Sen. *MicroRNA as biomarkers and diagnostics*. Journal of cellular physiology 231.1 (2016): 25-30.
- Hayes, Josie, Pier Paolo Peruzzi, and Sean Lawler. *MicroRNAs in cancer: biomarkers, functions and therapy*. Trends in molecular medicine 20.8 (2014): 460-469.
- Kappel, Andreas, et al. *MicroRNA in vitro diagnostics using immunoassay analyzers*. Clinical chemistry 61.4 (2015): 600-607.

miRNA quantification

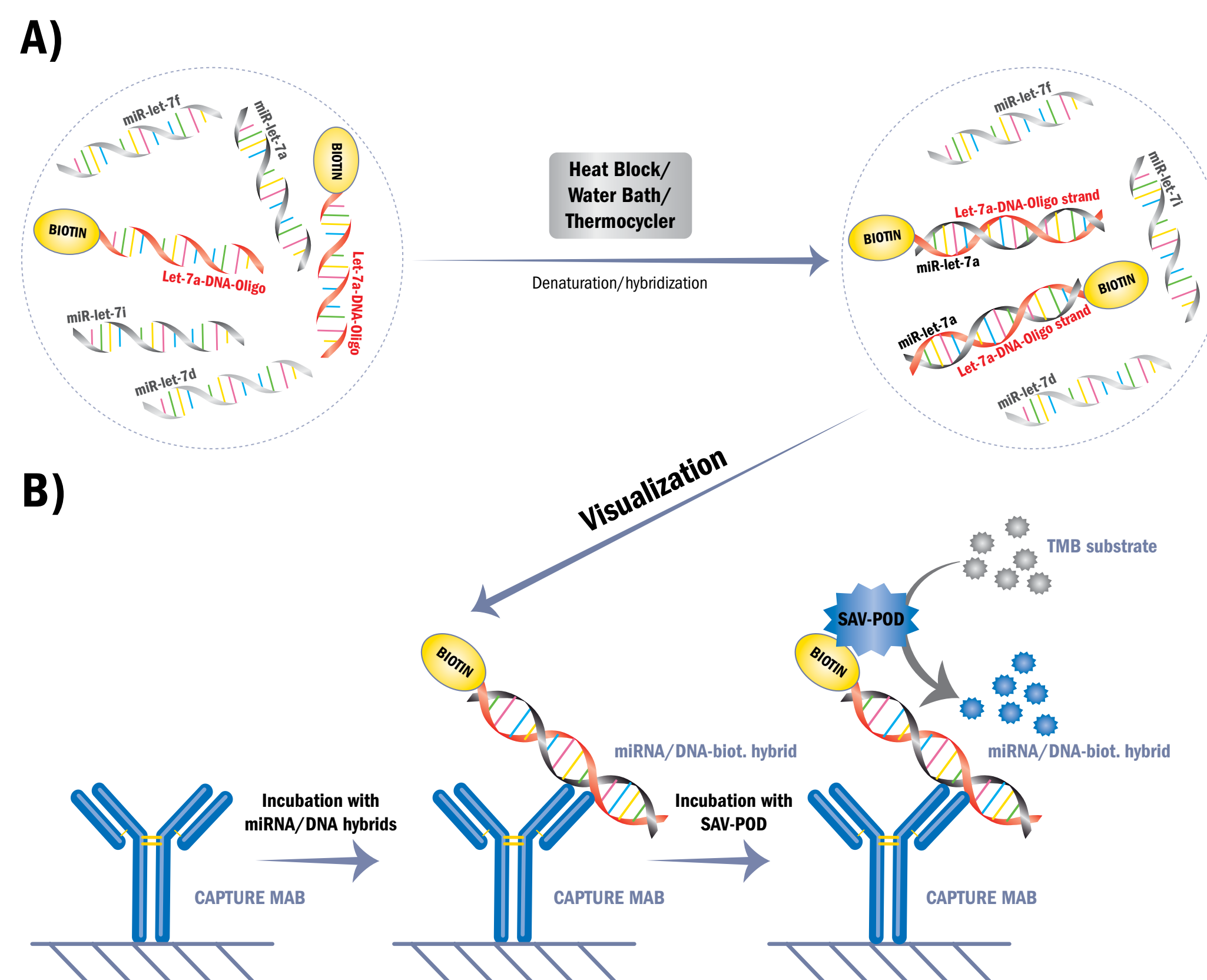

miREIA-21-5p


C) Calibration range

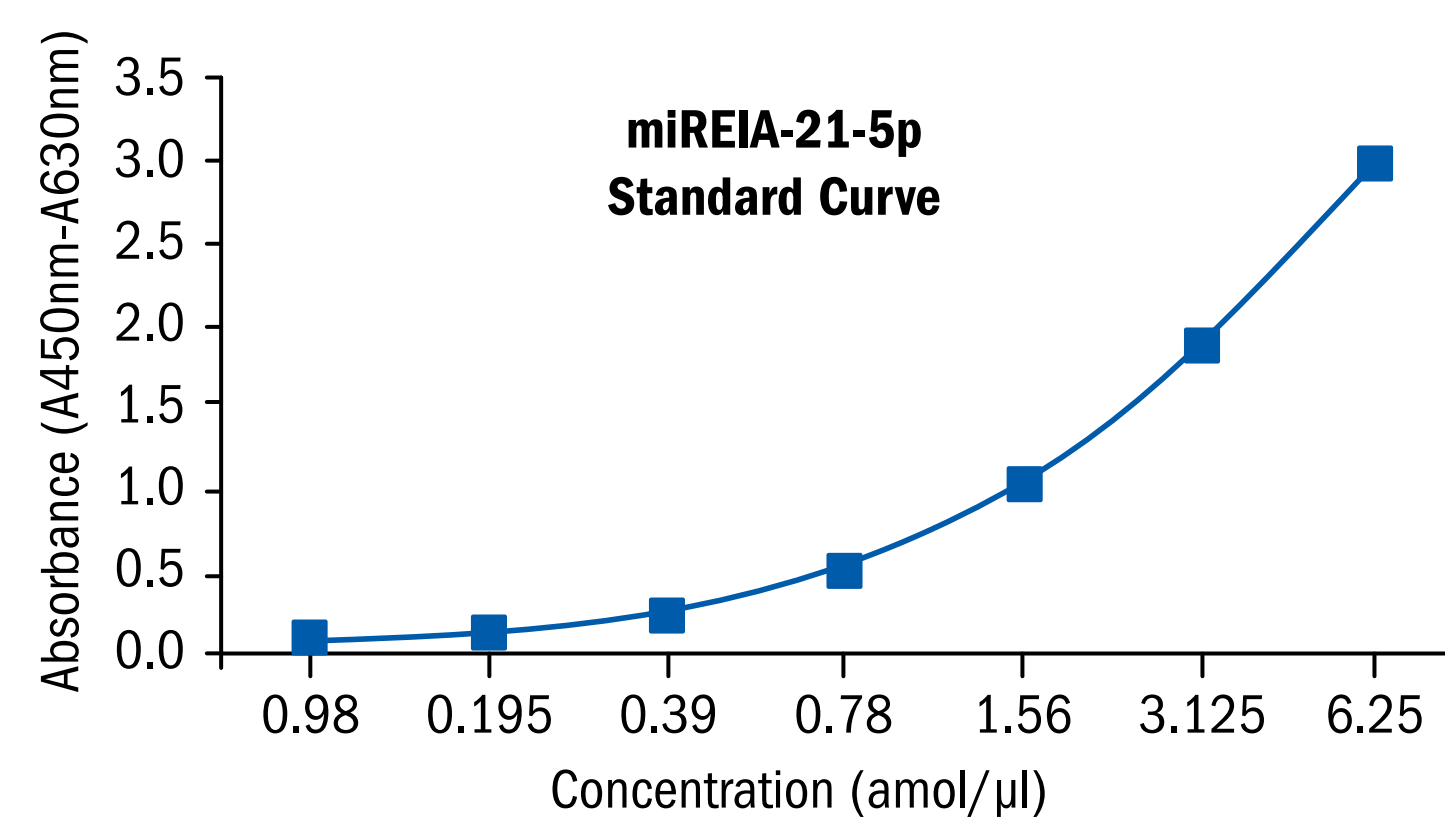

The calibration range of the miREIA-21-5p assay is 0.098-6.25 amol/ul.

E) Comparison with qRT-PCR

microRNA-21 isolated from whole blood taken from 10 healthy volunteers was quantified by miREIA-21-5p and by qRT-PCR using LNA nucleotide probes from Exiqon.

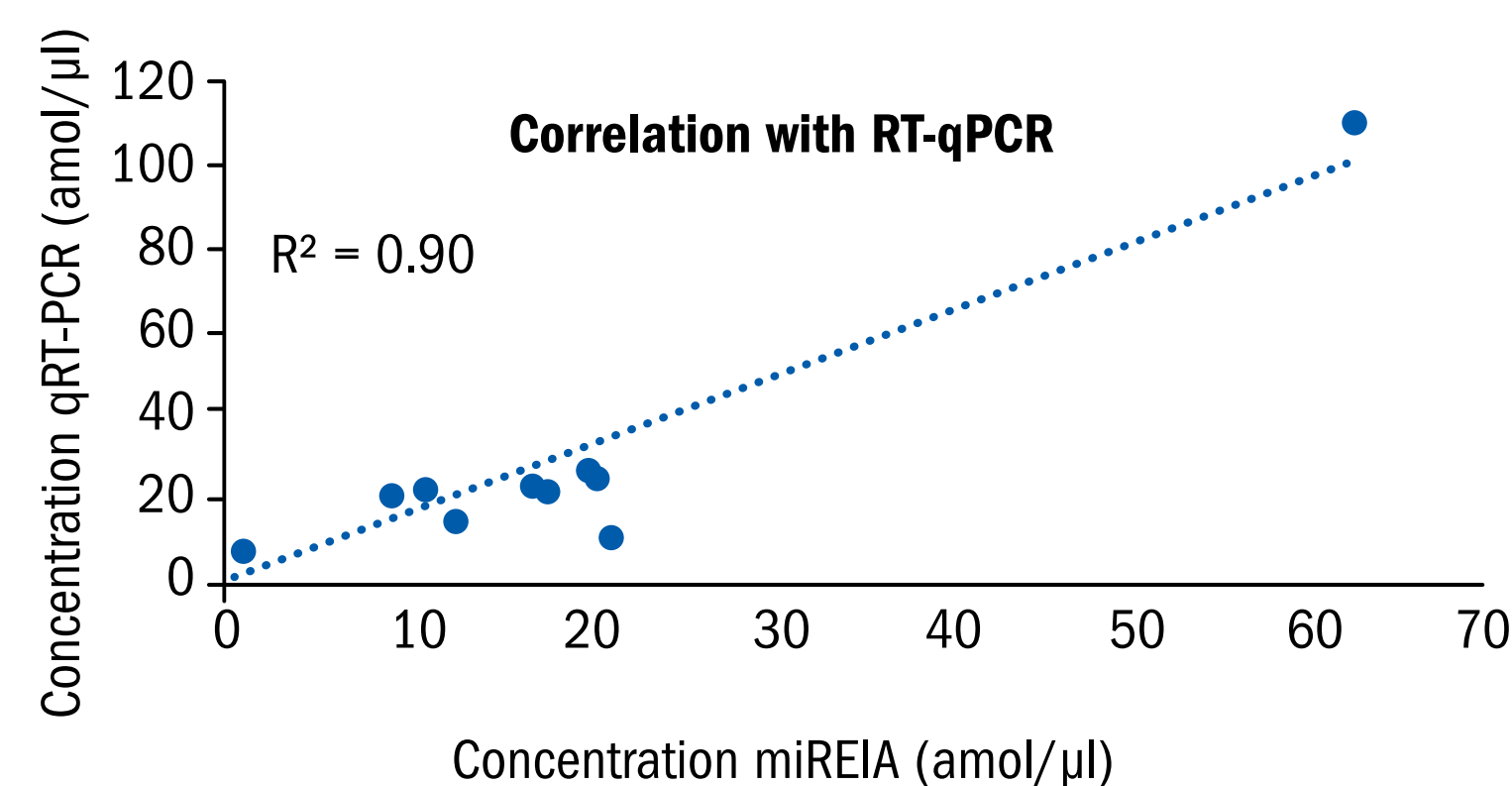

We found a strong correlation between qRT-PCR and miREIA-21-5p, the Pearson correlation coefficient being >0.9.


D) Dilution linearity

Two microRNA isolates from whole blood samples were serially diluted.

Sample	Sample dilution	Observed concentration amol/μl	Expected concentration amol/μl	Recovery O/E (%)
-	-	10.2	-	-
Sample1	2	5.0	5.1	99
	4	2.7	2.6	105
	8	1.3	1.3	100
-	-	20.4	-	-
Sample 2	2	10.1	10.2	99
	4	4.8	5.1	94
	8	2.6	2.6	100

Recovery was determined to be 94 - 105%

F) Total Assay Time

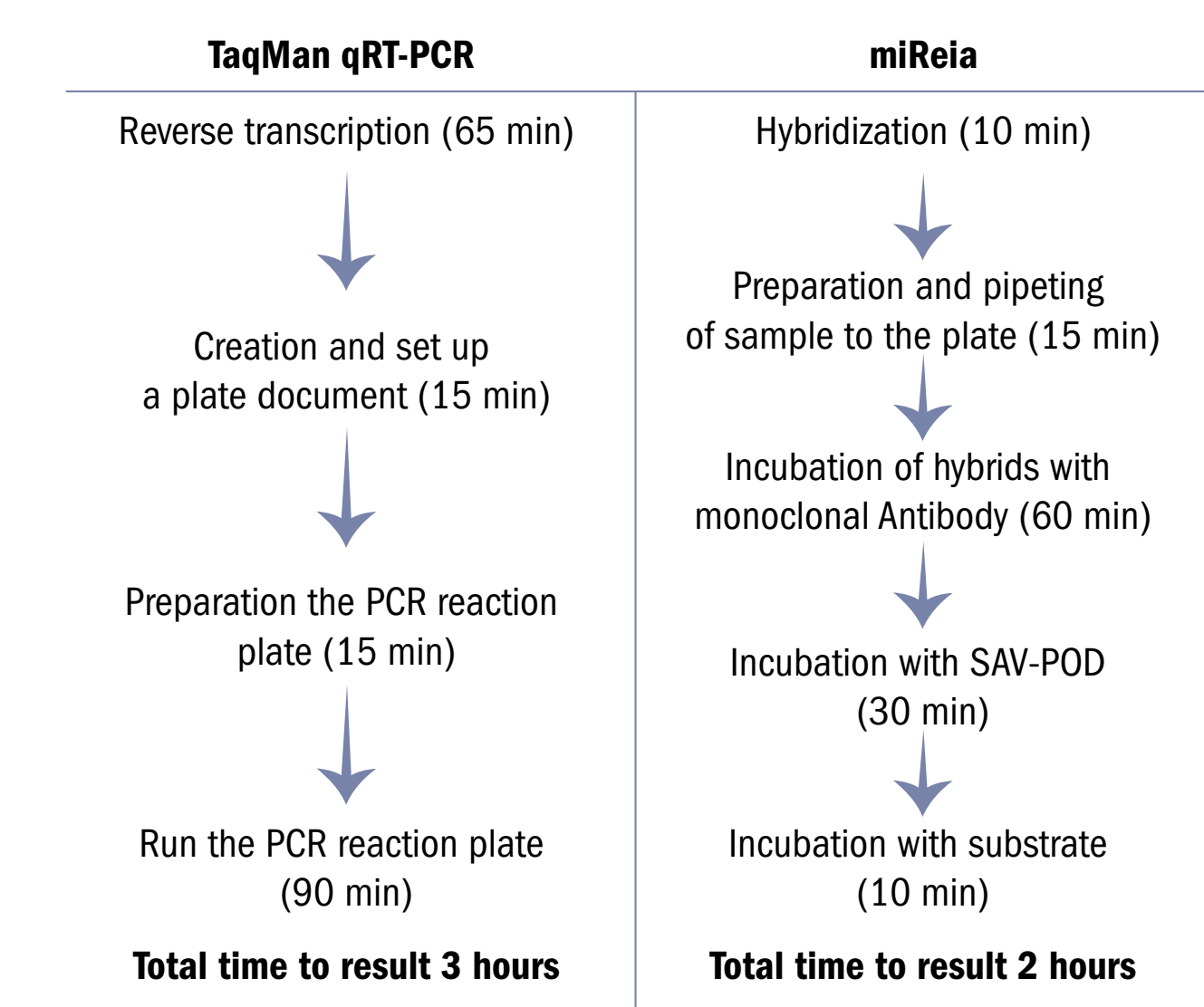

We compared the total time needed for miRNA quantification by two different methods: qRT-PCR and miREIA-21-5p. The total assay time of miREIA-21-5p is significantly shorter when compared to qRT-PCR.

Conclusion

- ▶ We are introducing a novel method for miRNA quantification
- ▶ miREIA-21-5p assay exhibits excellent analytical characteristics and strong correlation with qRT-PCR